


# Quick Guide

## FluoMini Pro Optical CF Seed Sensor

### Display & buttons

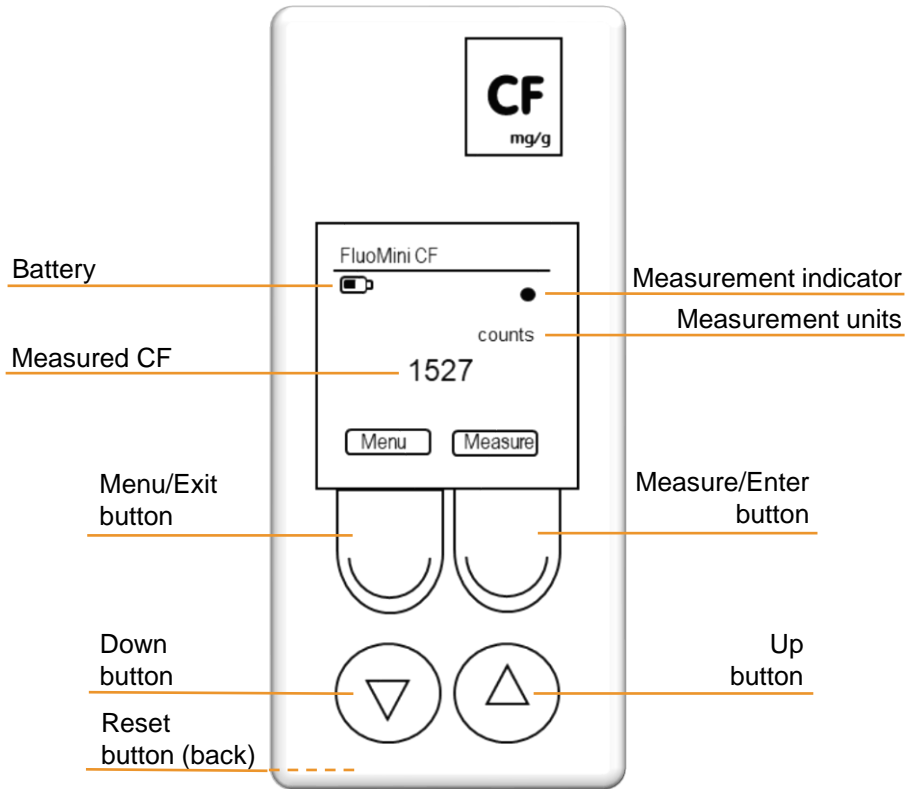

**Menu/Exit button**

Enter & exit main menu

**Measure/Enter button**

End sleep mode  
Single measurement > press briefly  
Start/stop continuous measurement > press 2 sec



# Sendot





www.sendot.nl

### Calibration

#### Apply reference sample


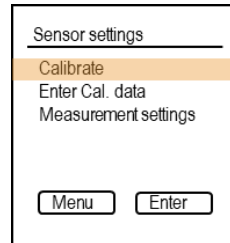
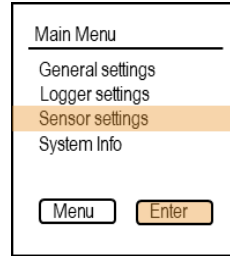


#### Choose between 1- or 2-point calibration


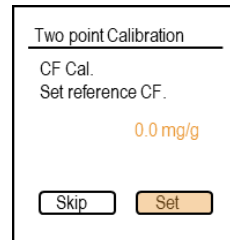
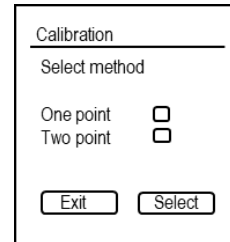


### Log function

#### Set interval time

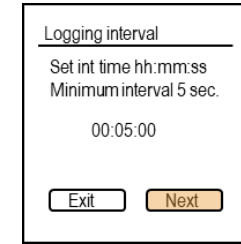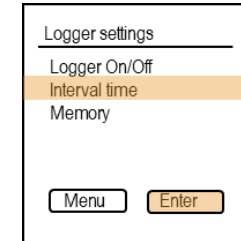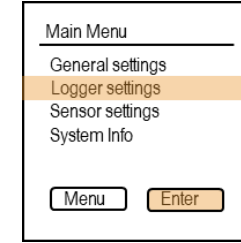



#### Start logging

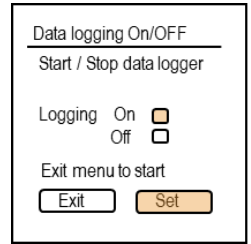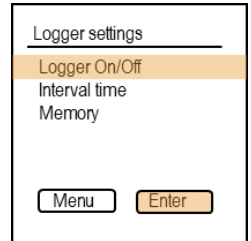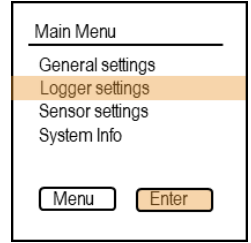

#### CAUTION!

Ensure to keep the glass of the probe holder clean to avoid falsified measurements. For further information read the manuals on our website.

#### Remove reference sample

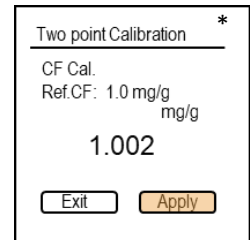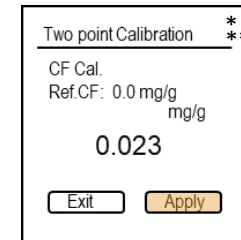

\* Wait for a stable measurement

\*\* Only for 2-point calibration